

## Neurokinin-1 (NK1) Human Receptor (Catalog #: RA25003) WB Protocol

- 1) Separate ~50 ug of whole cell or membrane protein on a 4-20% Tris-glycine gel.
- 2) Transfer the proteins from the gel onto a nitrocellulose membrane.
- 3) Incubate the nitrocellulose membrane in Blocking solution [5% non-fat dry milk / PBS] overnight at 4 degrees C.
- 4) Incubate the membrane with a 1:1,000 dilution of anti-NK-1 antibody (Novus Biologicals, Inc.,# NB300-119) in PBST-NFDM [PBS / 0.1% Tween-20 / 5% non-fat dry milk] for 2 hours at RT.
- 5) Wash the membrane with PBST-NFDM.
- 6) Incubate the membrane with HRP-conjugated goat anti-rabbit IgG diluted with PBST-NFDM for 1 hour at RT.
- 7) Detect antibody with an enhanced chemiluminescence reaction (ECL Western Blotting detection kit, Pharmacia).

## FACS Protocol

1. Seed cells into small (T-25) flask. Add 5-10 ml media. Allow cells to grow to 80-100% confluence.
2. When flasks are confluent, empty media and add new media containing appropriate drug, hormone, or cytokine dose.
3. Allow flasks to incubate for 24 hours. After incubation period, remove 2 ml of media and place 1 ml into one of two eppendorf tubes.
4. Empty off remaining media. Wash with 5 ml 1X PBS.
5. Add 2 ml PBS with azide (1% Goat serum). Scrape cells using a cell scraper.
6. Transfer cells to appropriate labeled FACSCAN tubes. Vortex all tubes
7. Incubate on ice for \_ hour in the PBS with azide (1% Goat serum). Vortex all tubes after incubation.
8. Centrifuge at 1400 RPM for 5 minutes. Empty off media.
9. Resuspend in 1 ml of a 1:3,000 primary antibody solution (NB 300-119) [1 ul antibody per 3 ml PBS with azide (1% Goat serum)].
10. Incubate on ice for 1 hour. Vortex all tubes after incubation.

---

FOR RESEARCH USE ONLY

Neuromics' reagents are for in vitro and certain non-human in vivo experimental use only and not intended for use in any human clinical investigation, diagnosis, prognosis, or treatment. We disclaim all liability in connection with the use of the information contained herein or otherwise, and all such risks are assumed by the user.

v1

11. Centrifuge at 1400 RPM for 5 minutes. Aspirate (or decant) supernatant.
12. Wash twice with 2 ml ice-cold PBS with azide (1% Goat serum).
13. Aspirate (or decant) supernatant and place tubes on ice.
14. Resuspend in 1 ml of a 1:500 secondary antibody solution [2 ul antibody per 1 ml PBS with azide (1% Goat serum)].
15. Vortex all tubes.
16. Cover with foil. Incubate for 30 minutes on ice.
17. Vortex after incubation. Centrifuge at 1400 RPM for 5 minutes.
18. Aspirate (or decant) supernatant. Wash twice with 2 ml ice-cold PBS with azide (1% Goat serum).
19. Aspirate (or decant) supernatant. Resuspend in 2% paraformaldehyde.
20. Read on flow.