



E18 Primary Dorsal Root Ganglia (DRGs)

Catalog#: PC35115

Components Included: 10 ganglion from one, embryonic day 1 from Sprague Dawley rat.

Storage Media-2 ml Hibernate A®/B27®/GlutaMAX™=**HAB**

Dissociation Media-5 ml Hibernate A without Calcium and Magnesium=**HA -Ca -Mg** **Culture and Growth Media-**12 ml NbActiv4 + NGF included.

Description: Primary Rat DRGs are live neurons isolated from micro-surgically dissected regions of day 18 embryonic Sprague/Dawley rat. These cells are prepared fresh each week and shipped in a nutrient rich medium that keeps the cells alive for up to 7 days under refrigeration. The cells are ideal for a wide variety of applications including: transfection, pharmacology, electrophysiology, immunocytochemistry, and neuronal development studies

Shipping/Storage: DRGs are shipped refrigerated. Cells are stable for up to 7 days when stored at 4-8°C. It is recommended to plate the cells as soon as possible after receiving cells according to Application Notes outlined below.

Protocol

Additional Components Required:

Dissociation Enzyme: Collagenase/Dispase (Roche 10 269 638 001).

Coating: Poly-D-Lysine (Sigma P6407).

Additional Growth Media: NbActiv4 + NGF Please contact Brett @ 866-350-1500 or brett@neuromics.com for additional media. You will need to add NGF (Invitrogen 13290-010) to the NbActiv4 (Catalog #: M36107).

Coverslips and dishes: 15mm round glass coverslips (Carolina Biological Supply #63-3031) and sterile 35mm petri dish (VWR# 25382-064).

Preparations (Room Temperature in a Sterile Hood):

1. Prepare substrate by coating with poly-D-lysine (100 µg/ml). Incubate coated surfaces for at least 1 hour (up to 20). Aspirate the poly-D-lysine, rinse once with ddH₂O, aspirate and air dry.
2. Prepare a 100 mg/ml stock solution of Collagenase/Dispase (Roche; Ref: 10 269 638 001) in ddH₂O.
3. To make the cell dissociation solution, dilute Step 2 to 1:100 with Hibernate A without Calcium and Magnesium (**HA -Ca -Mg**) for a final working solution of 1 mg/ml collagenase (0.1 U) / dispase (0.8 U). *Solution must be sterile filtered prior to use.* It can be aliquoted and frozen @ -20°C and thawed day of use.
4. Fire polish the tip of a sterile 9" silanized pasteur pipette to an opening of ~0.5 mm
5. Aliquot 20 µl of Trypan Blue (Sigma: T8154) into a 0.5 ml tube for Step 9.

Cell Dispersal (Room Temperature in a Sterile Hood):

1. With the silanized pasteur pipette, carefully remove the storage solution leaving the DRG's in minimal solution.
2. Place tissue in cell dissociation solution (Step 3, above) for 1 hr and incubate at 37°C (ambient atmosphere). Gently swirl every 5 min.
3. Allow tissue to settle and remove cell dissociation solution leaving the tissue at the bottom in minimal solution.
4. Immediately add 2 ml of Hibernate A®/B27®/GlutaMAX™ (**HAB**).
5. With the silanized pasteur pipette, triturate tissue ~30 times until 90% dispersed while avoiding air bubbles.
6. Let undispersed pieces settle for 1 min.
7. Transfer supernatant containing dispersed cells to a sterile 15 ml tube. Leave ~50 µl of HAB containing debris.

FOR RESEARCH USE ONLY V8-09/2014

Neuromics' reagents are for in vitro and certain non-human in vivo experimental use only and not intended for use in any human clinical investigation, diagnosis, prognosis, or treatment. We disclaim all liability in connection with the use of the information contained herein or otherwise, and all such risks are assumed by the user.

Neuromics • 5325 West 74th Street, Suite 8 • Edina, MN 55439 • phone 866-350-1500 • fax 612-677-3976 • e-mail pshuster@neuromics.com

8. Spin 1100 rpm (200 x G), 3 min. Discard supernatant leaving ~50 µl of HAB containing the pellet.
9. Disperse the pellet of cells (flick the bottom of the tube with a finger) and resuspend pellet in 1 ml NbActiv1® with 25 ng/mL NGF.
10. Aliquot 20 µl of cell solution into the 0.5 ml tube containing 20 µl of Trypan Blue (1:2 dilution).
11. Count cells using a hemacytometer (calculate cells/ml).

Cell Plating (Room Temperature in a Sterile Hood):

1. Dilute cells with NbActiv4® + 25 ng/ml NGF (0.2 ml/cm²) and plate at 100,000cells/cm² or desired concentration.
2. Incubate 37°C, 5% CO₂, 9% O₂, 95% humidity (or ambient O₂).
3. After 4 days, DRG's display axons and dendrites; synapses and action potentials begin at 7 days.
4. Change ½ of the medium with fresh, 37oC, CO2 equilibrated NbActiv4® + 25 ng/ml NGF every 3-4 days.

Viability Assay:

- Rinse twice with 37°C HBSS (0.2 ml/cm² of substrate).
- Prepare dye mix from an acetone stock of 15 mg/ml fluorescein diacetate and an aqueous stock of 4.6 mg/ml propidium iodide, dilute 15 µl of each into 1.5 ml HBSS (1:100 dilution).
- Add 20 µl of dye mix from step 2 to every 0.2 ml of HBSS added in step 1 (1:10 dilution).
- After ~1 min count live cells using blue excitation appropriate for fluorescein fluorescence (green cells). Count dead cells with green excitation for propidium iodide fluorescence (small red nuclei).
- Viability = (green cells/unit area)/(total cells plated/unit area) or Survival = green cells/(green + red cells)

References

A dissection and Tissue Culture Manual of the Nervous System (1989). A. Shahar, J.D. Vellis, A. Vernadakis, B. Haber (Eds.), *Dissociated Spinal Cord - Dorsal Root Ganglion Cultures on Plastic Tissue Culture Dishes and Glass Coverslips and Wells* (pp.219-222). Wiley-Liss, Inc.

J.L. Werth, S.A. Thayer (1994) Mitochondria Buffer Physiological Calcium Loads in Cultured Rat Dorsal Root Ganglion Neurons, *The Journal of Neuroscience*, 14(1), 348-356.

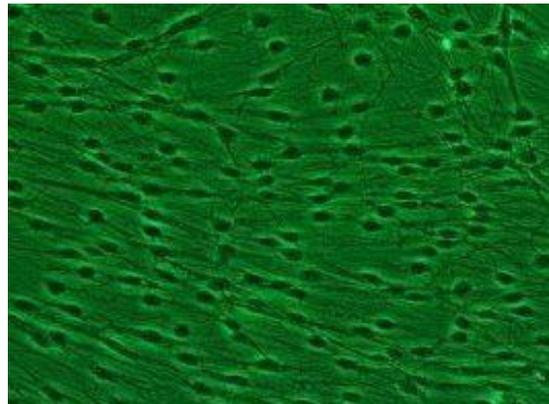


Image: DRGs on Poly-D-Lysine

FOR RESEARCH USE ONLY V8-09/2014

Neuromics' reagents are for in vitro and certain non-human in vivo experimental use only and not intended for use in any human clinical investigation, diagnosis, prognosis, or treatment. We disclaim all liability in connection with the use of the information contained herein or otherwise, and all such risks are assumed by the user.

Neuromics • 5325 West 74th Street, Suite 8 • Edina, MN 55439 • phone 866-350-1500 • fax 612-677-3976 • e-mail pshuster@neuromics.com