

# **Data Sheet**

# Gravity flow Strep-Tactin®XT Superflow® Column

Cat. No.: 2-4011-005, 2-4012-001,

2-4013-001, 2-4014-001

Lot No.:

Version: 6.0 Revision Date: 14.07.2020

IBA Headquarters IBA GmbH

Rudolf-Wissell-Str. 28 37079 Goettingen Germany

Tel. +49 (0) 551-5 06 72-0 Fax +49 (0) 551-5 06 72-181 IBA US contact information Fax 1-888-531-6813

E-mail: info@iba-lifesciences.com http://www.iba-lifesciences.com

| Description             | Immobilized streptavidin variant called <b>Strep-Tactin®XT</b> (5 mg/ml resin) which has been especially optimized for the purification of Strep-tag®II and Twin-Strep-tag® fusion proteins. |
|-------------------------|--|
| Support                 | Superflow® 6 (6% agarose, crosslinked)   |
| Form                    | Pre-packed in buffer:<br>100 mM Tris-HCl pH 8.0, 1.5 M (NH <sub>4</sub> ) <sub>2</sub> SO <sub>4</sub> , 150 mM NaCl,<br>1 mM EDTA, 0.02 % sodium azide.                                     |
| Biotin binding capacity | > 300 nmol/ml resin  |
| Stability               | 6 months after shipping  |
| Storage                 | recommended: 2-8 °C  |
| Shipping                | room temperature   |
| Hazards                 | Product is not classified as hazardous according to (EC) No 1272/2008 [CLP].  A Material Safety Data Sheet is provided.  |

| Elution      | Biotin Elution Buffer BXT (Buffer BXT):<br>100 mM Tris-HCl pH 8.0, 150 mM NaCl, 1 mM EDTA, 50 mM Biotin   |
|--------------|---|
| Regeneration | It is recommended to regenerate the column by using Strep-Tactin®XT Regeneration Buffer (3 M MgCl2, Buffer XT-R Cat. No. 2-1045-250). Alternatively, freshly prepared 10 mM NaOH can be used. |



Please download all up-to-date manuals, protocols and other material from <a href="https://www.iba-lifesciences.com/home.html">https://www.iba-lifesciences.com/home.html</a>.

## For research use only

### Important licensing information

This product is based on Strep-Tactin®XT technology covered by intellectual property (IP) rights and on completion of the sale IBA grants respective Limited Use Label Licenses to purchaser. IP rights and Limited Use Label Licenses for said technology are further described and identified at <a href="http://www.iba-lifesciences.com/patents.html">http://www.iba-lifesciences.com/patents.html</a> upon inquiry at <a href="mainto:info@iba-lifesciences.com">info@iba-lifesciences.com</a> or at IBA GmbH, Rudolf-Wissell-Str. 28, 37079 Goettingen, Germany. By use of this product the purchaser accepts the terms and conditions of all applicable Limited Use Label Licenses.

### Trademark information