

# Data Sheet

## StrepMan Magnet

### for 15ml and 50ml tubes

Cat. No.: 6-5650-065

Last date of revision  
March 2020  
Version 5650-1

<b>Description</b>	<b>StrepMan permanent magnet for 2x15 ml and 2x50 ml conical reaction tubes. To be used for manual cell isolation with Strep-Tactin® Magnetic Microbeads using the Streptamer® Technology.</b>
<b>Attention</b>	<p><b>Strong permanent magnet.</b></p> <p><b>The Strep-Tactin® Magnetic Microbeads (MB) have to be stored distant from strong magnetic fields like the StrepMan Magnet (at least 10 cm).</b></p> <p><b>Danger of injury by strong magnetic attraction of ferromagnetic materials such as iron.</b></p> <ul style="list-style-type: none"> <li>• <b>Keep away from magnetic storage devices such as hard disks, diskettes, ID cards, credit cards, otherwise stored data may be deleted.</b></li> <li>• <b>Keep away from electronic devices and persons with cardiac pacemakers.</b></li> <li>• <b>Persons with cardiac pacemakers should not be in close contact with this magnet.</b></li> </ul>
<b>Cave!</b>	<b>For better handling, a lubricant should be applied to the rubber rings from time to time. Then the tubes can be better inserted and the rubber rings do not go off.</b>
<b>Cleaning</b>	<b>Use only commercially available cleaning agents such as mild detergent or alcohol (70% ethanol) to clean the StrepMan. Do not autoclave.</b>
<b>Dimensions</b>	<b>W 13,6 × D 9,7 × H 12 cm; Weight 900 g</b>
<b>Warranty</b>	<b>24 months</b>



Go digital and help the environment. Please download all up-to-date manuals, protocols and other material from <http://www.iba-lifesciences.com>.

#### For research use only

##### Important licensing information

This product is based on Twin-Strep-tag® and Streptamer® technologies covered by intellectual property (IP) rights and on completion of the sale IBA grants respective Limited Use Label Licenses to purchaser. IP rights and Limited Use Label Licenses for said technology are further described and identified at <http://www.iba-lifesciences.com/patents.html> or upon inquiry at [info@iba-lifesciences.com](mailto:info@iba-lifesciences.com) or at IBA GmbH, Rudolf-Wissell-Str. 28, 37079 Goettingen, Germany. By use of this product the purchaser accepts the terms and conditions of all applicable Limited Use Label Licenses.

##### Trademark information

The owners of trademarks marked by “®” or “TM” are identified at <http://www.iba-lifesciences.com/patents.html>. Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.